

INSTRUCTION SHEET


Transpector[®] MPH/MPS Installation Instructions 961-416-P1/P2 Heating Jacket

Introduction

The Transpector MPH/MPS Heater Jacket Mount Kit (961-433-G1) provided with Heating Jacket Kit (961-029) is intended to inhibit the movement of the Manifold Heater. This kit is only for system installations without calibration standards that are mounted in a vertical orientation.

NOTE: For all other heater installations, discard this kit.

- 1** Install the Manifold Heater over the tube extension and align the snaps with the tapped hole in the sensor flange.
- 2** Align the Internal Heater Bracket so the large hole fits over the snap on the heating jacket and capture the bracket in the Manifold Heater by securing the snaps.
- 3** Using a 7 mm combination wrench, attach the Internal Heater Bracket to the tapped hole in the sensor flange using the hex head fastener provided in the kit.
- 4** Attach the manifold heater power connector to the mating connector on the 600-1117 power cable provided in the 961-029 heating jacket kit.

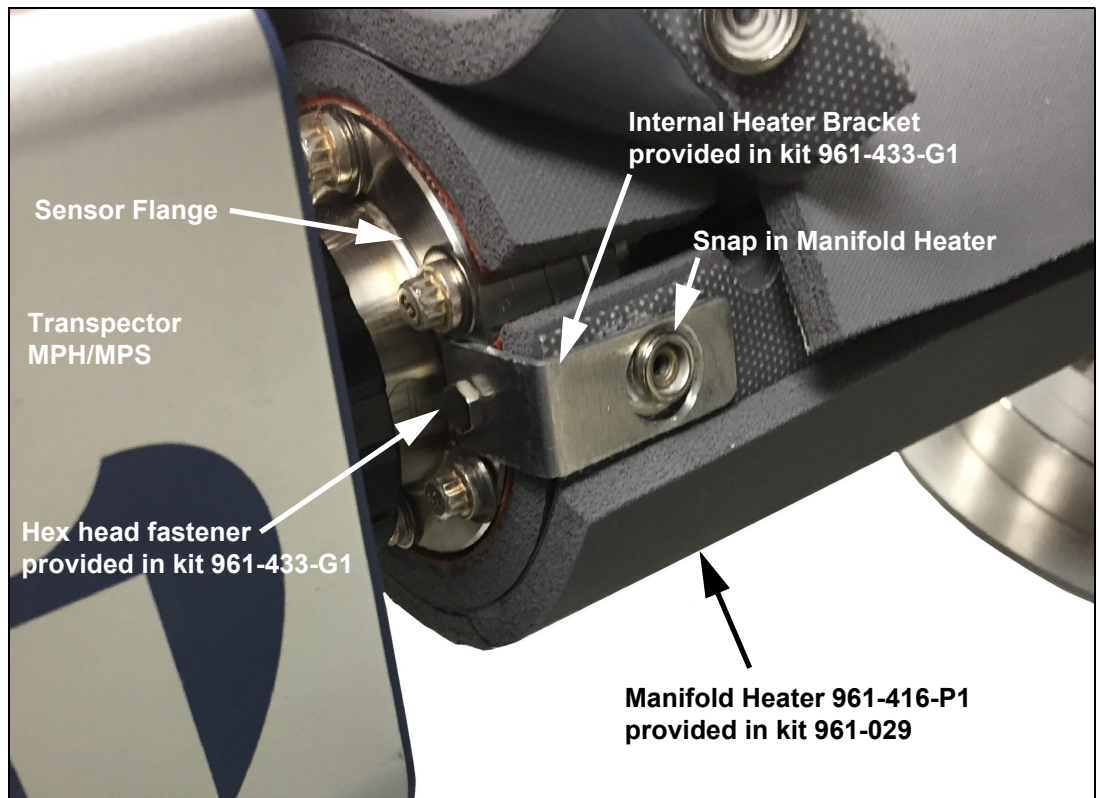

www.inficon.com reachus@inficon.com

Due to our continuing program of product improvements, specifications are subject to change without notice. All trademarks are the property of their respective owners.

IPN 074-679-P1A ©2016 INFICON